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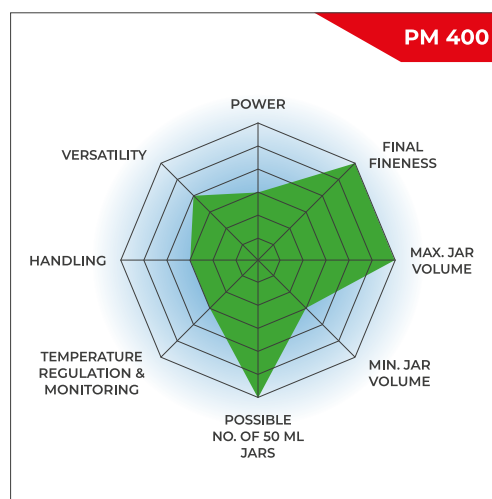
THE HIGH-VOLUME BALL MILL FOR HIGH THROUGHPUT APPLICATIONS

- | Max. speed 400 rpm, large sun wheel
- | Up to 10 mm feed size and 0.1 µm final fineness
- | 4 grinding stations for jars from 12 ml up to 500 ml, jars of 12 – 80 ml can be stacked (two jars each)
- | GrindControl to measure temperature and pressure inside the jar.
- | Aeration lids to control the atmosphere inside the jar
- | Floor model, storable SOPs and cycle programs, 5 different jar materials for dry and wet grinding

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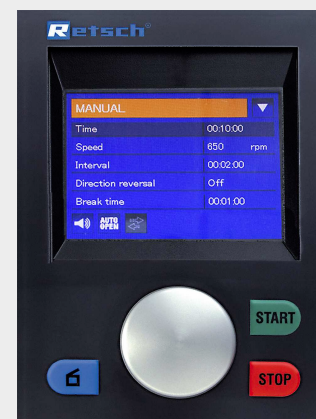
THE BEST ALTERNATIVE TO A RETSCH PLANETARY BALL MILL? A RETSCH MIXER MILL.

Benefit from particularly ergonomic handling while achieving the same finenesses down to the nanometer range.

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EASYFIT GRINDING JARS FOR EXCELLENT RESULTS

The performance and the result of sample preparation are also determined by the choice of the grinding jar and its ball charge. The EasyFit range of jars has been specially designed for extreme working conditions such as long-term trials, even at maximum speed of 800 rpm, wet grinding, high mechanical loads and maximum speeds as well as for mechanical alloying. This line of jars is suitable for all RETSCH planetary ball mills.

The new EasyFit grinding jar series features a structure on the bottom of the 50-500 ml jars called Advanced Anti-Twist (AAT). This ensures that the jars are tightly fixed without the risk of twisting, even at high speed, and that wear and tear is drastically reduced. Secure clamping of the jars is made much easier: to find the correct clamping position, a maximum twist of 60° is required.

The geometry of the EasyFit jars in the 50 ml and 250 ml sizes has been enlarged in diameter and reduced in height compared to the previous "comfort" models. This offers two advantages: better grinding results and interchangeable lids, as there are only three diameter dimensions for the entire grinding jar range.

Diameter categories

- | Diameter 1: 12 ml and 25 ml grinding jars
- | Diameter 2: 50 ml, 80 ml and 125 ml grinding jars
- | Diameter 3: 250 ml and 500 ml grinding jars

- | Available jar sizes: 12 ml / 25 ml / 50 ml / 80 ml / 125 ml / 250 ml / 500 ml
- | Innovative Advanced Anti-Twist (AAT) function ensures secure fit of grinding jars
- | High flexibility thanks to suitability of three lid sizes for all seven jar sizes
- | Pressure-tight and dust-proof O-ring sealing prevents material spillage
- | Jars and balls available in 5 materials: hardened stainless steel, tungsten carbide, agate, sintered aluminium oxide, zirconium oxide
- | Stainless steel protective jacket for agate, sintered aluminum oxide, zirconium oxide and tungsten carbide grinding jars
- | Зазор между стаканом и крышкой позволяет легко открыть крышку, например, с помощью шпателя, если внутри стакана возникает пониженное давление

JARS & LIDS FOR SPECIAL APPLICATIONS

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Both the aeration lid and GrindControl can now be equipped with inlays of different materials. Thus, the lid can be used for, e. g. a steel and a zirconium oxide jar by simply exchanging the inlay.

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Video:
Aeration lid

ADAPTER FOR SPECIAL APPLICATIONS

With a special adapter, co-crystal screening can be carried out in a planetary ball mill, using disposable vials such as 1.5 ml GC glass vials. The adapter features 24 positions arranged in an outer ring with 16 positions and an inner ring with 8 positions. The outer ring accepts up to 16 vials, allowing for screening up to 64 samples simultaneously when using the Planetary Ball Mill PM 400. The 8 positions of the inner ring are suitable to perform trials with different energy input, e.g. for mechanosynthesis research.

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22.661.0002  Clamping unit for PM 100 / PM 400

03.025.0178 Adapter for stacking grinding jars 50 ml - 80 ml

02.728.0048  Counter aid for sun wheel PM 100, PM 200 and PM 400

03.486.0062 Opening aid for clamping unit of planetary ball mills

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PRESSURE AND TEMPERATURE MEASURING SYSTEM GRINDCONTROL FOR PLANETARY BALL MILLS

**incl. sensors and transmitter unit, insert of lid, software, case, opening aid and cleaning accessories for PM
(please order grinding jars separately)**

22.782.0033

GrindControl for PM grinding jar EasyFit 50 - 125 ml

22.782.0034

GrindControl for PM grinding jar EasyFit 250 - 500 ml

GRINDCONTROL LID INSERTS

03.474.0243 GrindControl lid insert for 50, 80, 125 ml, stainless steel

03.474.0246 GrindControl lid insert for 50, 80, 125 ml, zirconium oxide

03.474.0244 GrindControl lid insert for 250 or 500 ml, stainless steel

03.474.0247 GrindControl lid insert for 250 or 500 ml, zirconium oxide

ACCESSORIES FOR PM GRINDCONTROL WITH GRINDING JARS EASYFIT

05.114.0056

O-ring for 50, 80 or 125 ml

05.114.0054

O-ring for 250 ml - 500 ml grinding jars EasyFit (PM)

03.111.0438 Flat gasket for 50 ml, 80 ml or 125 ml

03.111.0439 Flat gasket for 250 ml - 500 ml

22.186.0007 Sintered filter with O-ring, set of 10 pieces

22.864.0001

Valve set M8x1 for GrindControl and aeration lids

GRINDING JARS EASYFIT

(grinding jars EasyFit are suitable for all planetary ball mills)

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ADAPTER FOR GLASS VIALS

01.462.0540		Adapter for 24 x 1.5 ml glass vials, stainless, hardened steel
22.749.0009		Glass vial 1.5 ml incl. septum cap, 100 pieces
05.181.0112		Replacement pressure spring for adapter for 24 x 1.5 ml glass vials, 1 piece
01.462.0541		Adapter for 7 x 20 ml glass vials, stainless, hardened steel
22.749.0010		Glass vial 20 ml incl. septum cap, 100 pieces
05.181.0044		Replacement pressure spring for adapter for 7 x 20 ml glass vials, 1 piece

ACCESSORIES FOR GRINDING JARS EASYFIT FOR WET GRINDING, GRINDING WITH INERT ATMOSPHERE AND MECHANICAL ALLOYING (MA)

AERATION LIDS (INCL. INLAY)

22.107.0613	for grinding jars EasyFit 50 ml - 125 ml, hardened stainless steel
22.107.0616	for grinding jars EasyFit 50 ml - 125 ml, tungsten carbide
22.107.0617	for grinding jars EasyFit 50 ml - 125 ml, agate
22.107.0615	for grinding jars EasyFit 50 ml - 125 ml, zirconium oxide
22.107.0618	for grinding jars EasyFit 250 ml - 500 ml, hardened stainless steel
22.107.0621	for grinding jars EasyFit 250 ml - 500 ml, tungsten carbide
22.107.0622	for grinding jars EasyFit 250 ml - 500 ml, agate
22.107.0620	for grinding jars EasyFit 250 ml - 500 ml, zirconium oxide
22.107.0619	for grinding jars EasyFit 250 ml - 500 ml, aluminum oxide
22.864.0001	 Spare valve set for aeration lids M8x1

INLAY FOR AERATION LID

03.474.0225	for grinding jars EasyFit 50 ml - 125 ml, hardened stainless steel
03.474.0207	for grinding jars EasyFit 50 ml - 125 ml, tungsten carbide
03.474.0208	for grinding jars EasyFit 50 ml - 125 ml, agate
03.474.0206	for grinding jars EasyFit 50 ml - 125 ml, zirconium oxide
03.474.0226	for grinding jars EasyFit 250 ml - 500 ml, hardened stainless steel
03.474.0210	for grinding jars EasyFit 250 ml - 500 ml, tungsten carbide
03.474.0211	for grinding jars EasyFit 250 ml - 500 ml, agate
03.474.0209	for grinding jars EasyFit 250 ml - 500 ml, zirconium oxide
03.474.0215	for grinding jars EasyFit 250 ml - 500 ml, aluminum oxide

AERATION LIDS FOR GRINDING JARS EASYFIT

INCL. O-RINGS AND SINTERED FILTER (PLEASE ORDER LID INSERT AND GRINDING JAR SEPARATELY)

22.107.0636	Aeration lid for grinding jar EasyFit 50 ml - 125 ml
22.107.0637	Aeration lid for grinding jar EasyFit 250 ml - 500 ml

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22.867.0011	for grinding jars EasyFit 50 ml - 125 ml
22.867.0012	for grinding jars EasyFit 250 ml - 500 ml
02.486.0055	Opening aid for safety closure device

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