XRD-MOLEN MCCRONE

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ADVANTAGES

- | crystal lattice preserved during grinding operation
- | very narrow and reproducible particle size distribution
- | minimal cross contamination
- | compact, bench-top sized model
- | adjustable grinding power (4 steps)
- | geschikt voor droog en nat malen
- | easy to clean
- | almost maintenance-free
- | quiet operation

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beenderen, boriden, bouwmaterialen, carbiden, cement, glas, keramiek, klei, leisteen, metalen, mica, mineralen, nitriden, plantenmaterialen

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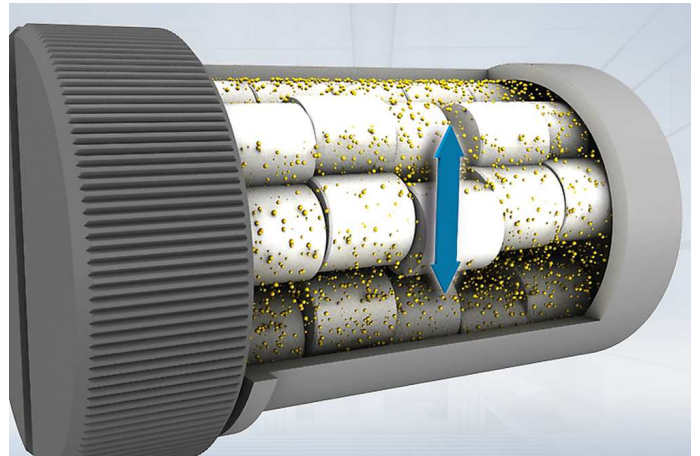
PRINCIPE

The XRD-Mill McCrone carries out size reduction mainly by friction. 48 cylindrical grinding elements are placed into the grinding jar in eight rows of six elements each.

During operation, the jar's circular motion causes the elements to grind the sample from < 0.5 mm to the low μm -range (typically < 10 μm).

Thanks to the very gentle size reduction process, the crystal lattice of the sample is preserved. This makes the XRD-Mill McCrone the instrument of choice for the sample preparation for subsequent X-ray diffraction analysis.

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BESTELGEGEVENS

XRD-MILL MCCRONE

**XRD-Mill McCrone with 2 grinding jars with screw top lid of PP, 1 pouring lid, 1 loading device
(Please order grinding elements separately)**

20.770.0001

XRD-Mill McCrone, 220-240 V, 50/60 Hz

XRD-MILL MCCRONE COMPLETE PACKAGE

XRD-Mill McCrone complete package incl. 2 agate grinding sets (50 grinding elements each), 2 grinding jars with screw top lid of PP, 1 pouring lid, 1 loading device

20.770.1001

XRD-Mill McCrone, 220-240 V, 50/60 Hz complete package

SETS OF GRINDING ELEMENTS

02.368.0123

Agaat

1 set (50 grinding elements)

02.368.0122

Corundum

1 set (48 grinding elements)

02.368.0126

Zirconium oxide, 1 set (48 grinding elements)

ACCESSORIES XRD-MILL MCCRONE

22.460.0001

Sample preparation set (stainless steel percussion mortar, 10 corundum grinding elements, 1 sieve 500 µm and 1 brush)

02.368.0124

Grinding elements, tungsten carbide, 2 pieces

ADDITIONAL ITEMS XRD-MILL MCCRONE

01.462.0356

Grinding jar with screw top lid

03.462.0359

Screw top lid

03.462.0357

Pouring lid

03.486.0007

Loading device

22.111.0005

Gasket for grinding jar, 5 pcs.